




Application Note
Pharmaceutical
USP Class 1 & 2
Residual Solvents



www.defiant-tech.com

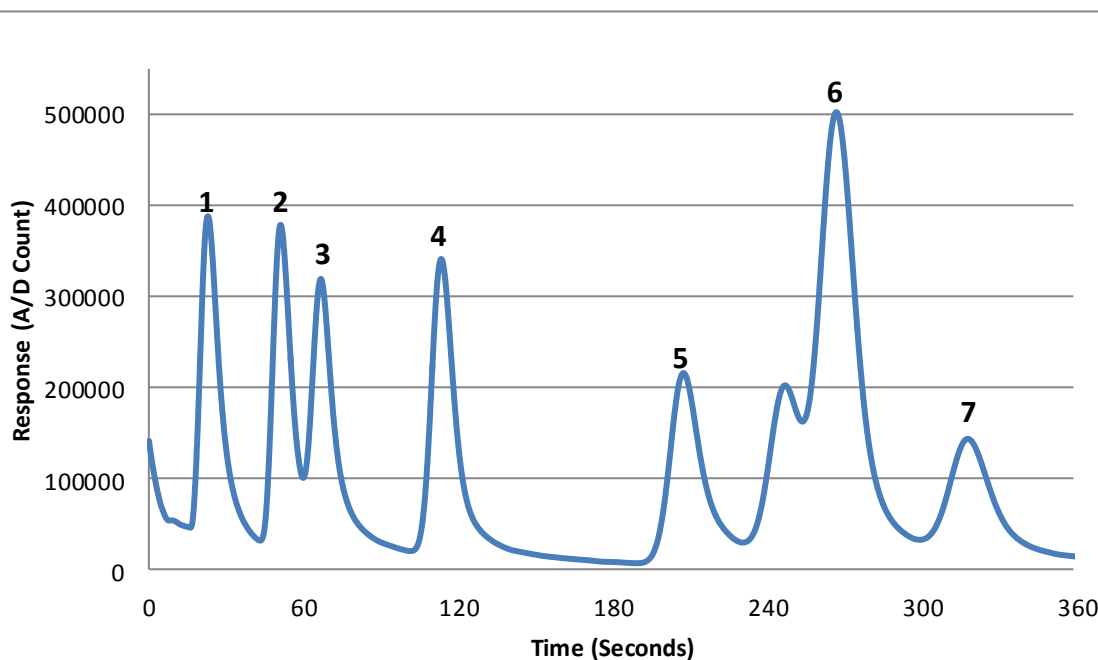

Hand held micro system for the **6-minute** detection of:

- 1) 1,1-Dichloroethene
- 2) Benzene
- 3) Trichloroethylene
- 4) Toluene
- 5) Chlorobenzene
- 6) p/m-Xylene
- 7) o-Xylene

FROG-4000™

Gas Chromatography System for VOCs

Instrument Settings:

2.5m Column, 0.80 Micron OV-1 Deposition

Collect 30 seconds

GC Ramp

Cold 40°C hold for 360 seconds

Ramp 40°C to 60°C over 60 seconds

Hot 100°C hold for 30 seconds

Instrument Specifications:

Weight: 4.8lbs.

Dimensions: 10X7.5X14.5inches

Battery Life: 6 Hours, or External Power

Interface: USB Port or Bluetooth



Defiant Technologies Inc.
6814A Academy Parkway W NE
Albuquerque, NM 87109
505-999-5880



www.defiant-tech.com