




Application Note
Pharmaceutical
USP Class 3 Residual
Solvents




www.defiant-tech.com

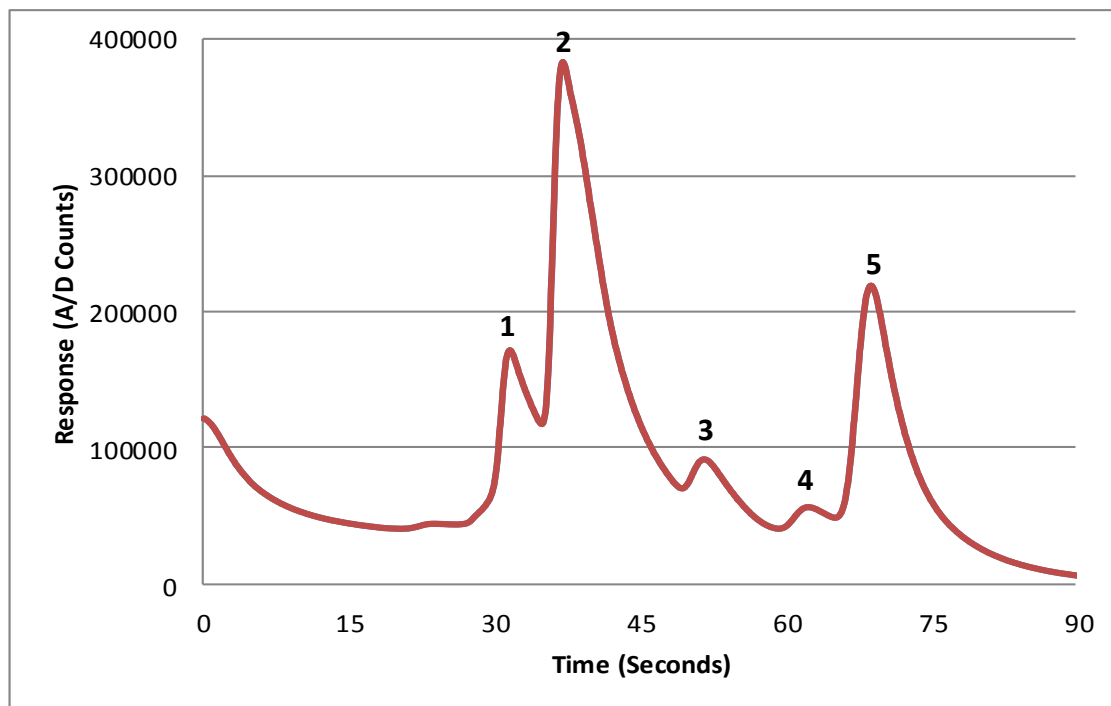

Hand held micro system for the detection of:

- 1) Acetone
- 2) Ethanol
- 3) Ethyl Acetate
- 4) 1-Propanol
- 5) 2-Propanol

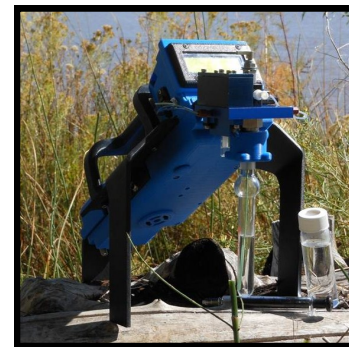

FROG-4000™

Gas Chromatography System for VOCs

Instrument Settings:

2.5m Column, 0.80 Micron OV-1 Deposition

Collect 30 seconds

GC Ramp

Cold 40°C hold for 90 seconds

Ramp 40°C to 60°C over 30 seconds

Hot 100°C hold for 60 seconds

Instrument Specifications:

Weight: 4.8lbs.

Dimensions: 10X7.5X14.5inches

Battery Life: 6 Hours, or External Power

Interface: USB Port or Bluetooth



Defiant Technologies Inc.
6814A Academy Parkway W NE
Albuquerque, NM 87109
505-999-5880



www.defiant-tech.com