



FROG-4000
Installing the Sparge
Bottle Ferrule



This is the correct installation for the sparge bottle ferrule. Correct installation is crucial for proper operation.



The pictures on the left and right show the proper installation of the sparge bottle ferrule. On the left the ferrule is shown on a globe neck sparge bottle. On the right, it is shown installed on an air sample adapter.



If your retention times aren't lining up correctly or you cannot get a consistent calibration. First inspect the sparge bottle ferrule. If it is improperly installed, it can create a small leak in the system and cause this variation.